Centricon® Centrifugal Filter Device

— Product Overview —

For complete information about this product, please contact Millipore Corporation.

Usage Guidelines

- Centricon devices with Ultracel® YM-10, -30, -50, and -100 membranes are approved for *in-vitro* diagnostic use.
- Centricon devices with Ultracel YM-3 membranes are for research use only; not for use in diagnostic procedures.
- Do not autoclave Centricon components.
- Do not operate above the following limits.

Centricon-3 devices	without retentate vial:	6500 × g
	with retentate vial:	7500 × g
	during prerinsing:	5000 × g
Centricon-10, -30, and -50 devices:		5000 × g
Centricon-100 devices	with proteins:	5000 × g
	with nucleic acids:	2000 × g
- The ultrafiltration membranes in Centricon devices contain trace amounts of glycerine (~2 µL per device). Prerinse if necessary. Do not allow the membrane to dry out once wet.

Instructions for Use

After inserting sample reservoir into filtrate vial, add solution to sample reservoir. Place covered device into centrifuge and counterbalance with similar device. Spin until desired concentration is achieved. To recover retentate, place retentate vial over sample reservoir, invert unit, and centrifuge at 300–1000 × g for 2 minutes.

Specifications

Sample reservoir: styrene acrylonitrile
Membrane support base: polycarbonate
O-Ring: red silicone rubber
Filtrate and retentate vials, cap: polypropylene
Membrane: Ultracel YM regenerated cellulose membrane
Maximum initial sample volume: 2.0 mL
Typical final concentration volume: 25–50 µL
Hold-up volume: 0.02 mL
Active membrane area: 0.92 cm²

Technical Assistance

Find up-to-date worldwide contact information at www.millipore.com/offices.

Warranty

This product is covered by the standard limited warranty, as set forth in Millipore's Life Science Catalogue.

P36573, Rev. —, 08/04

© 2004, Millipore Corporation. All rights reserved. Printed in U.S.A.

Millipore, Centricon, and Ultracel are registered trademarks of Millipore Corporation.

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